



# UNPACKING INSTRUCTIONS

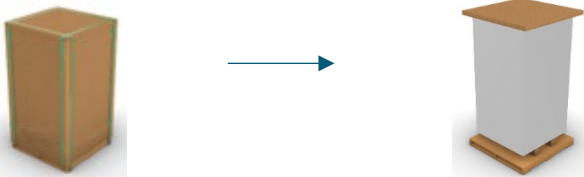
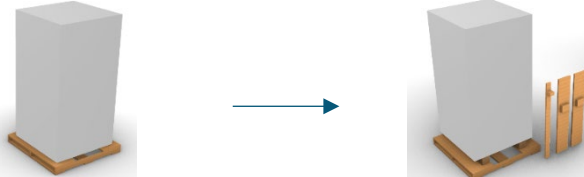

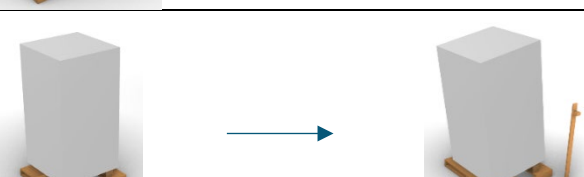


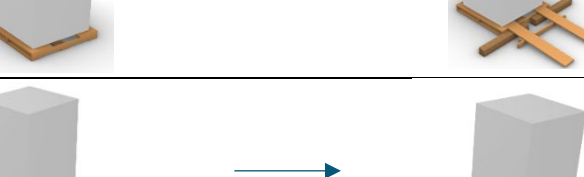
Large Upright Freezers



Scan for Video



## RECOMMENDED TOOLS: Scissors & Phillips Screwdriver

Step	Illustration	Instructions
1		Remove straps, plastic screws, top cardboard-frame, cardboard sides, foam, and utility bag from unit.
2		Remove wood lever and ramps from within pallet. Ensure all casters are oriented in the same direction and parallel to guide rail. Stow ramps somewhere accessible for step 6.
3		Use lever to prop-up unit.
4		While unit is propped, remove one corrugated cardboard buildup from underneath unit. Remove lever. Save the buildup for step 6.
5		Repeat steps 3 & 4 to remove the remaining corrugated cardboard buildup from underneath unit.
6		Set up ramps on side of pallet by lining up the brackets on the ramps with the notches on the pallet. Use both corrugated cardboard buildups and lever to add support to the ramps.
7		Carefully maneuver unit down the ramps using three to four people.

Upon receiving the unit, we advise you to properly dispose of any packaging material used to secure the unit during shipping. Please refer to local recycling codes and dispose of materials accordingly. For information on waste disposal sites, please refer to your local waste management company or recycling center.

Global Cooling technology is manufactured under U.S. and International patents. Stirling Ultracold is a trademark of Global Cooling, Inc.

**STIRLING POWER**

[www.stirlingultracold.com](http://www.stirlingultracold.com)